

Saccharomyces cerevisiae Host Cell Proteins

Immunoenzymetric Assay for the Measurement of Saccharomyces cerevisiae Host Cell Proteins Catalog # F135

Intended Use

This kit is intended for use in determining the presence of *Saccharomyces cerevisiae* protein impurities in products manufactured by recombinant expression in this yeast cell line. The kit is for **Research and Manufacturing Use Only** and is not intended for diagnostic use in humans or animals.

Summary and Explanation

Recombinant expression of proteins by *S.cerevisiae* is an efficient method to obtain cost effective quantities of a desired protein. Many of these recombinantly produced proteins are intended for use as therapeutic agents in humans and animals and as such must be highly purified. The manufacturing and purification process of these products leaves the potential for impurities by host cell proteins from *S.cerevisiae*. Such impurities can result in adverse toxic or immunological reactions and thus it is desirable to reduce host cell impurity to the lowest levels practical.

Immunological methods using antibodies to HCPs such as Western Blot and ELISA are conventionally accepted. While Western blot is a useful method aiding in the identity of HCPs, it suffers from a number of limitations. Western blot is a complex and technique dependent procedure requiring a subjective interpretation of results. Furthermore, it is essentially a qualitative method and does not lend itself to obtaining quantitative answers. The sensitivity of Western blot is severely limited by the volume of sample that can be tested and by interference from the presence of high concentrations of the intended product. Western Blot may be able to detect HCPs in samples from upstream in the purification process but it often lacks adequate sensitivity and specificity to detect HCPs in purified downstream and final product. The microtiter plate immunoenzymetric assay (ELISA) method employed in this kit overcomes the limitations of Western blots providing on the order of 100 fold better sensitivity. This simple to use, highly sensitive, objective, and semi-quantitative ELISA is a powerful method to aid in optimal purification process development, process control, routine quality control, and product release testing. This kit is "generic" in the sense that it is intended to react with essentially all of the HCPs that could contaminate the product independent of the purification process. The

antibodies have been generated against and affinity purified using a mild lysate washed of *S.cerevisiae* cells to obtain HCPs typically encountered in your initial product recovery step. Western blot was used as a preliminary method and established that the antibodies reacted to the majority of HCP bands resolved by the PAGE separation. If you have need of a more sensitive and specific method to demonstrate reactivity to individual HCPs in your samples *Cygnus Technologies* recommends a method we find superior to 2D Western blot. If you have a need of a more sensitive method to demonstrate coverage to HCPs in your process *Cygnus Technologies* recommends a method that is superior to Western blot called Antibody Affinity Extraction (AAE). AAE is has greatly increased sensitivity and specificity to Western blot which makes it a better predictor of how the antibodies will perform in the ELISA. For additional information on AAE please visit our website and read the posted articles under Technical Documents or contact our Technical Services Department.

Special procedures were utilized in the generation of these antibodies to ensure that low molecular weight and less immunogenic impurities as well as high molecular weight components would be represented. As such this kit can be used as a process development tool to monitor the optimal removal of host cell impurities as well as in routine final product release testing. Because of the high sensitivity and broad reactivity of the antibodies, this generic kit has been successfully qualified for testing of final product HCPs in many different products regardless of growth and purification process. When the kit can be satisfactorily qualified for your samples, the application of a more process specific assay is probably not necessary in that such an assay would only provide information redundant to this generic assay. However, if your qualification studies indicate the antibodies in this kit are not sufficiently reactive with your process specific HCPs it may be desirable to also develop a more process specific ELISA. This later generation assay may require the use of a more specific and defined antisera. Alternatively, if the polyclonal antibody used in this kit provides sufficient sensitivity and broad antigen reactivity, it may be possible to substitute the standards used in this kit for ones made from the impurities that typically co-purify through your purification process and thus achieve better accuracy for process specific HCPs. The suitability of this kit for a given sample type and product must be determined and

qualified experimentally by each laboratory. The use of a process specific assay with more defined antigens and antibodies in theory may yield better sensitivity however such an assay runs the risk of being too specific in that it may fail to detect new or atypical impurities that might result from some process irregularity or change. For this reason it is recommended that a broadly reactive "generic" host cell protein assay be used as part of the final product purity analysis even when a process specific assay is available. If you deem a more process specific assay is necessary, *Cygnus Technologies* is available to apply its proven technologies to develop such antibodies and assays on custom basis.

Reagents & Materials Provided

Component	Product #
Anti-<i>S.cerevisiae</i>, biotinylated Affinity purified goat antibody conjugated to biotin in a protein matrix with preservative. 1x12mL	F136
Anti-<i>S.cerevisiae</i> coated microtiter strips 12x8 well strips in a bag with desiccant	F137*
<i>S.cerevisiae</i> Standards Solubilized <i>S.cerevisiae</i> HCPs in bovine serum albumin with preservative. Standards at 0, 2, 8, 25, 75, and 200ng/mL. 1 mL/vial	F138
Streptavidin:Alkaline Phosphatase In a protein matrix with preservative. 1x12mL.	F009B
PNPP Substrate p-nitrophenyl phosphate in a Diethanolamine buffer with preservative. 1x12mL.	F008
Wash Concentrate (20X) Tris buffered saline with preservative. 1x50mL.	F004

*All components can be purchased separately except # F137.

Principle of the Procedure

The *Saccharomyces cerevisiae* Host Cell Protein assay is a two-site immunoenzymetric assay. Samples containing *S.cerevisiae* proteins are reacted in microtiter strips coated with an affinity purified anti-*S.cerevisiae* capture antibody. A second biotinylated anti-*S.cerevisiae* antibody is reacted simultaneously forming a sandwich complex of solid phase antibody-*S.cerevisiae* protein-biotin labeled antibody. The microtiter strips are then washed to remove any unbound reactants. Streptavidin:Alkaline Phosphatase is added to bind to any biotinylated antibody bound to the well. The substrate p-nitrophenyl phosphate (PNPP) is then reacted. The amount of hydrolyzed substrate is read on a microtiter plate reader and will be directly proportional to the concentration of *S.cerevisiae* proteins present.

Storage & Stability

- All reagents should be stored at 2°C to 8°C for stability until the expiration date printed on the kit.
- The substrate reagent should not be used if its absorbance at 405nm is greater than 0.4.
- Reconstituted wash solution is stable until the expiration date of the kit.
- After prolonged storage, you may notice a salt precipitate and/or yellowing of the wash concentrate. These changes will not impact assay performance. To dissolve the precipitate, mix the wash concentrate thoroughly and dilute as directed in the 'Preparation of Reagents' section.

Materials & Equipment Required But Not Provided

- Microtiter plate reader spectrophotometer with dual wavelength capability at 405 & 492nm. (*If your plate reader does not provide dual wavelength analysis you may read at just the 405nm wavelength.*)
- Pipettors - 50µL and 100µL
- Repeating or multichannel pipettor - 100µL
- Microtiter plate rotator (400 - 600 rpm)
- Sample Diluent (recommended Cat # 1028)
- Distilled water
- 1 liter wash bottle for diluted wash solution

Precautions

- For Research or Manufacturing use only.
- This kit should only be used by qualified technicians.

Preparation of Reagents

- Bring all reagents to room temperature.
- Dilute wash concentrate to 1 liter in distilled water, label with kit lot and expiration date, and store at 4°C.

Procedural Notes

1. Complete washing of the plates to remove excess unreacted reagents is essential to good assay reproducibility and sensitivity. We advise against the use of automated or other manual operated vacuum aspiration devices for washing plates as these may result in lower specific absorbances, higher non-specific absorbance, and more variable precision. The manual wash procedure described below generally provides lower backgrounds, higher specific absorbance, and

better precision. If duplicate CVs are poor or if the absorbance of the 0 standard is greater than 0.300, evaluate plate washing procedure for proper performance.

2. High Dose Hook Effect or poor dilutional linearity may be observed in samples with very high concentrations of HCP. High Dose Hook Effect is due to insufficient excess of antibody for very high concentrations of HCPs present in the samples upstream in the purification process. Samples greater than 50 µg/mL may give absorbances less than the 200 ng/mL standard. It is also possible for samples to have certain HCPs in concentrations exceeding the amount of antibody for that particular HCP. In such cases the absorbance of the undiluted sample may be lower than the highest standard in the kit however these samples will fail to show acceptable dilutional linearity/parallelism as evidenced by an apparent increase in dilution corrected HCP concentration with increasing dilution. High Dose Hook and poor dilutional linearity are most likely to be encountered from samples early in the purification process. If a hook effect is possible, samples should also be assayed diluted. If the HCP concentration of the undiluted sample is less than the diluted sample this may be indicative of the hook effect. Such samples should be diluted at least to the minimum required dilutions (MRDs) as established by your qualification studies using your actual final and in-process drug samples. The MRD is the first dilution at which all subsequent dilutions yield the same HCP value within the statistical limits of assay precision. The HCP value to be reported for such samples is the dilution corrected value at or greater than the established MRD. The diluent used should be compatible with accurate recovery. The preferred diluent is our Cat# 1028 available in 100mL, 500mL, or 1 liter bottles. This is the same material used to prepare the kit standards. As the sample is diluted in 1028 its matrix begins to approach that of the standards thus reducing any inaccuracies caused by dilutional artifacts. Other prospective diluents must be tested for recovery by using them to dilute the 200ng/mL standard, as described in the "Limitations" section below.

3. If the substrate has a distinct yellow color prior to performing the assay it may have been contaminated. If this appears to be the case read 200µL of substrate against a water blank. If the absorbance is greater than 0.4 it may be necessary to obtain new substrate or the sensitivity of the assay may be compromised. The PNPP substrate is very sensitive to environmental impurities. Do not leave bottle open or at room temperature for longer than is needed. Only remove as much reagent as is needed for your assay run and do not return any unused substrate back into the substrate bottle. Additional substrate can be purchased separately as Cat # F008.

- The assay protocol is a simultaneous incubation of sample with biotinylated antibody. This yields a sensitivity of 0.5ng/mL and requires 4 hours to complete.
- The assay is very robust such that assay variables like incubation times, sample size, and other sequential incubation schemes can be altered to manipulate assay performance for more sensitivity, increased upper analytical range, or reduced sample matrix interference. Before modifying the protocol from what is recommended, users are advised to contact our technical services for input on the best way to achieve your desired goals.
- The protocol specifies the use of an approved orbital microtiter plate shaker for the immunological step. These can be purchased from most laboratory supply companies. If you do not have such a device it is possible to incubate the plate without shaking, however it will be necessary to extend the first immunological incubation step by one hour in order to achieve comparable results to the 2-hour shaking protocol. **Do not shake during the 1-hour substrate incubation step as this may result in higher backgrounds and worse precision.**
- Bring all reagents to room temperature.
- Set-up plate spectrophotometer to read dual wavelength at 405nm for the test wavelength and 492nm for the reference wavelength.
- All standards, controls and test samples should be assayed at least in duplicate.
- Maintain a repetitive timing sequence from well to well for all assay steps to ensure that all incubation times are the same for each well.
- Make a work list for each assay to identify the location of each standard control and patient.
- Thorough washing is essential to proper performance of this assay. Automated plate washing systems or other vacuum aspiration devices are not recommended. The manual method described in the assay protocol is preferred for best precision, sensitivity and accuracy. A more detailed discussion of this procedure can be obtained from our Technical Services Department or on our web site. In addition, a video demonstration of proper plate washing technique is available in the 'Technical Help' section of our web site.

Assay Protocol

1. Pipette 50 μ L of standards, controls and samples into wells indicated on work list.
2. Pipette 100 μ L of biotinylated anti-*S.cerevisiae* (#F136) into each well.
3. Cover & incubate on orbital shaker at 400 – 600 rpm for 2 hours at room temperature, 24°C \pm 4°C.
4. Dump contents of wells into waste. Blot and gently but firmly tap over absorbent paper to remove most of the residual liquid. Overly aggressive banging of the plate or use of vacuum aspiration devices in an attempt to remove all residual liquid is not necessary and may cause variable dissociation of antibody bound material resulting in lower ODs and worse precision. Fill wells generously to overflowing with diluted wash solution using a squirt bottle or by pipetting in ~350 μ L. Dump and tap again. Repeat for a total of 4 washes. Wipe off any liquid from the bottom outside of the microtiter wells as any residue can interfere in the reading step. Do not allow wash solution to remain in wells for longer than a few seconds. Do not allow wells to dry before adding substrate.
5. Pipette 100 μ L of Streptavidin: Alkaline Phosphatase (#F009B) into each well.
6. Cover & incubate on orbital shaker at 400 – 600 rpm for 1 hour at room temperature, 24°C \pm 4°C.
7. Repeat Step 4 as shown above.
8. Pipette 100 μ L of PNPP substrate (#F008).
9. Incubate for 1 hour at room temperature.
10. Read absorbance at 405/492nm.

Limitations

- Before relying exclusively on this assay to detect host cell proteins, each laboratory should qualify that the kit antibodies and assay procedure yield acceptable specificity, accuracy, and precision. A suggested protocol for this qualification can be obtained by contacting our Technical Services Department or at our web site.
- The standards used in this assay are comprised of *S.cerevisiae* HCPs solubilized by mechanical disruption and detergent. 1D Western blot analysis of the antibodies used in this kit demonstrates that they recognize the majority of distinct PAGE separated bands seen using a sensitive protein staining method like silver stain or colloidal gold. Because the vast majority of HCPs will be conserved among all strains of *S.cerevisiae*, this kit should be adequately reactive to HCPs from your strain. Other clients have successfully qualified this kit for their individual

S.cerevisiae strains demonstrating acceptable specificity, accuracy, and sensitivity for process intermediate samples as well as final product. However, there can be no guarantee that this assay will detect all proteins or protein fragments from your process. If you desire a much more sensitive and specific method than western blot to detect the reactivity of the antibodies in this kit to your individual HCPs *Cygnus* is pleased to perform AAE as a service to provide coverage information of the antibodies to the HCPs in your process samples.

- Certain sample matrices may interfere in this assay. However, the potential exists that the product protein itself, high or low pH, high salt, detergents, or other components in the sample matrix may result in either positive or negative interference in this assay. It is advised to test all sample matrices for interference by diluting the 200ng/mL standard, 1 part to 4 parts of the matrix containing no or very low HCP impurities. This diluted standard when assayed as an unknown should give a value of 30 to 50 ng/mL. Consult *Cygnus Technologies* Technical Service Department for advice on how to quantitate the assay in problematic matrices.

Calculation of Results

The standards may be used to construct a standard curve with values reported in ng/mL “total immunoreactive HCP equivalents” (See Limitations section above). This data reduction may be performed through computer methods using curve fitting routines such as point-to-point, cubic spline, or 4 parameter logistic fit. **Do not use linear regression analysis to interpolate values for samples as this may lead to significant inaccuracies!** Data may also be manually reduced by plotting the absorbance values of the standard on the y-axis versus concentration on the x-axis and drawing a smooth point-to-point line. Absorbances of samples are then interpolated from this standard curve.

Quality Control

- Precision on duplicate samples should yield average % coefficients of variation of less than 10% for samples greater than 2 ng/mL and less than 200 ng/mL. CVs for samples less than 2 ng/mL may be greater than 10%.
- For optimal performance the absorbance of the substrate when blanked against water should be less than 0.4.
- It is recommended that each laboratory assay appropriate quality control samples in each run to ensure that all reagents and procedures are correct. **You are strongly urged to make controls in your typical sample matrix using HCPs derived from**

your cell line. These controls can be aliquoted into single use vials and stored frozen for long-term stability.

variation (CVs) are less than 20%. The LOQ is 0.9ng/mL.

Precision

Both intra (n=20 replicates) and inter-assay (n=5 assays) precision were determined on 3 pools with low (2ng/mL), medium (75ng/mL), and high concentrations (200ng/mL). The % CV is the standard deviation divided by the mean and multiplied by 100.

Pool	Intra assay CV	Inter assay CV
Low	15.4%	16.1%
Medium	4.2%	5.6%
High	2.1%	3.4%

Specificity/Cross-Reactivity

1D Western blot analysis against other strains of *S.cerevisiae* indicates that most of the proteins are conserved among all strains. Thus, this assay should be useful for detecting HCP's from other *Saccharomyces cerevisiae* strains. Each end user must qualify that this kit is adequately reactive and specific for their samples. 1D Western blot is highly orthogonal to ELISA and to non-specific protein staining methods such as silver stain or colloidal gold. As such, the lack of identity between silver stain and western blot does not necessarily mean there is no antibody to that protein or that the ELISA will not detect that protein. If you desire a much more sensitive and specific method than Western blot to detect the reactivity of the antibodies in this kit to your individual HCPs *Cygnus* is pleased to perform AAE as a service to provide coverage information of the antibodies to the HCPs in your process samples. This method has been shown to be much at least 100 fold more sensitive than Western blots in detecting antibody reactivity to individual HCPs. The same antibody as is used for both capture and HRP label can be purchased separately.

Cell proteins from the yeast *Pichia pastoris* were tested and found to be non-reactive.

Recovery/ Interference Studies

Various buffer matrices were evaluated by adding known amounts of *S.cerevisiae* HCP preparation used to make the standards in this kit. Because this assay is designed to minimize matrix interference most of these buffers yielded acceptable recovery defined as between 80-120%. In general extremes in pH (less than 5.0 and greater than 8.5) as well as certain detergents can cause under-recovery. High concentrations of the recombinant product protein may also cause positive or negative interference in this assay. Each user should qualify that their sample matrices yield accurate recovery. Such an experiment can be performed, by diluting the 200ng/mL standard provided with this kit, into the sample matrix in question and assaying for recovery.

Example Data

Well #	Contents	Abs. at 405-490nm	Mean Abs.
A1	Zero Std	0.000	0.002
B1	Zero Std	0.004	
C1	2ng/mL	0.025	0.024
D1	2ng/mL	0.023	
E1	8ng/mL	0.101	0.098
F1	8ng/mL	0.095	
G1	25ng/mL	0.306	0.300
H1	25ng/mL	0.295	
A2	75ng/mL	0.879	0.872
B2	75ng/mL	0.865	
C2	200ng/mL	1.702	1.712
D2	200ng/mL	1.722	

Performance Characteristics

Cygnus Technologies has qualified this assay by conventional criteria as indicated below. A more detailed copy of this "Qualification Summary" report can be obtained by request. This qualification is generic in nature and is intended to supplement but not replace certain user and product specific qualification and qualification which should be performed by each laboratory. At a minimum each laboratory is urged to perform a spike and recovery study in their sample types. In addition, any of your samples types containing process derived HCPs within or above the analytical range of this assay should be evaluated for dilutional linearity to ensure that the assay is accurate and has sufficient antibody excess for your particular HCPs. Each laboratory and technician should also demonstrate competency in the assay by performing a precision study similar to that described below. A more detailed discussion of recommended user qualification protocols can be obtained by contacting our Technical Services Department or at our web site.

Sensitivity

The lower limit of detection (**LOD**) is defined as that concentration corresponding to a signal two standard deviations above the mean of the zero standard. LOD is 0.5ng/mL.

The lower limit of quantitation (**LOQ**) is defined as the lowest concentration, where concentration coefficients of

Hook Capacity

Increasing concentrations of HCPs greater than 200 ng/mL were assayed as unknowns. The hook capacity, defined as that concentration which will give an absorbance reading less than the 200 ng/mL standard was greater than 50 µg/mL.

Ordering Information/ Customer Service

Cygnus Technologies also offers kits for the extraction and detection of CHO Host Cell DNA. The following kits are available:

- Residual Host Cell DNA extraction:
Cat # D100W, DNA Extraction Kit in 96 deep well plate
Cat # D100T, DNA Extraction Kit in microfuge tubes

To place an order or to obtain additional product information contact *Cygnus Technologies*:

www.cygnustechnologies.com

Cygnus Technologies, LLC
4332 Southport Supply Rd. SE
Southport, NC 28461 USA
Tel: 910-454-9442
Fax: 910-454-9443
Email: techsupport@cygnustechnologies.com
