

E. coli Host Cell Proteins

Immunoenzymetric Assay for the Measurement of *E. coli* Host Cell Proteins Catalog # F410

Intended Use

This kit is intended for use in determining the presence of *E. coli* host cell protein impurities in products manufactured by recombinant expression in *E. coli*. The kit is for **Research and Manufacturing Use Only** and is not intended for diagnostic use in humans or animals.

Summary and Explanation

Recombinant expression in *E. coli* is a relatively simple and cost effective method for production of certain proteins and pDNA. Many of these products are intended for use as therapeutic agents in humans and animals, and as such, must be highly purified. The manufacturing and purification process of these products leaves the potential for impurities by host cell proteins (HCPs) from *E. coli*. Such impurities can reduce the efficacy of the therapeutic agent and result in adverse toxic or immunological reactions and thus it is desirable to reduce HCP impurities to the lowest levels practical.

Immunological methods using antibodies to HCPs such as Western Blot and ELISA are conventionally accepted. While Western blot is a useful method aiding in the identity of HCPs, it suffers from a number of limitations. Western blot is a complex and technique dependent procedure requiring subjective interpretation of results. Furthermore, it is essentially a qualitative method and does not lend itself to obtaining quantitative answers. The sensitivity of Western blot is severely limited by the volume of sample that can be tested and by interference from the presence of high concentrations of the intended product. While Western Blot may be able to detect HCPs in samples from upstream in the purification process, it often lacks adequate sensitivity and specificity to detect HCPs in purified downstream and final product. The microtiter plate immunoenzymetric assay (ELISA) method employed in this kit overcomes the limitations of Western blots providing on the order of 100 fold better sensitivity. This simple to use, highly sensitive, objective, and semi-quantitative ELISA is a powerful method to aid in optimal purification process development, process control, routine quality control, and product release testing. This kit is "generic" in the sense that it is intended to react with essentially all of the HCPs that could pollute the product independent of the purification process. The antibodies have been generated against and affinity purified using a blended lysate of E. coli from a variety of strains including the

commonly used DH5a and BL21 strains. This relatively mild lysing procedure is intended to obtain HCPs typically encountered in initial product recovery steps, such as clarification of conditioned media when the product is secreted, or after osmotic shock or mild detergent and mechanical disruption, to obtain inclusion bodies and other intracellular proteins. Western blot was used as a preliminary method and established that the antibodies reacted to the majority of HCP bands resolved by the PAGE separation. If you have a need of a more sensitive method to demonstrate coverage to HCPs in your process Cyanus Technologies recommends a method that is superior to Western blot called Antibody Affinity Extraction (AAE). AAE has greatly increased sensitivity and specificity to Western blot which makes it a better predictor of how the antibodies will perform in the ELISA. For additional information on AAE please visit our website and read the posted articles under Technical Documents or contact our Technical Services Department.

Special procedures were utilized in the generation of these antibodies to ensure that low molecular weight and less immunogenic impurities as well as high molecular weight components would be represented. As such, this kit can be used as a process development tool to monitor the optimal removal of host cell impurities as well as in routine final product release testing.

Because of the high sensitivity and broad reactivity of the antibodies, this generic kit has been successfully qualified for testing of final product HCPs in many different products regardless of growth and purification process. When the kit can be satisfactorily qualified for your samples, the application of a more process specific assay may not be necessary, in that such an assay would only provide information redundant to this generic assay. However, if your qualification studies indicate the antibodies in this kit are not sufficiently reactive with your process specific HCPs it may be desirable to also develop a more process specific ELISA. This later generation assay may require the use of a more specific and defined antisera. Alternatively, if the polyclonal antibody used in this kit provides sufficient sensitivity and broad antigen reactivity, it may be possible to substitute the standards used in this kit for ones made from the impurities that typically co-purify through your purification process and thus achieve better accuracy for process specific HCPs. The suitability of this kit for a given sample type and product must be determined and qualified experimentally by each laboratory. The use of a process specific assay with more defined antigens and antibodies in theory may yield better sensitivity however such an assay runs the risk of being too specific in that it may fail to detect new or atypical impurities that might result from some process irregularity or change. For this reason it is recommended that a broadly reactive "generic" host cell protein assay be used as part of the final product purity analysis even when a process specific assay is available. If you deem a more process specific assay is necessary, Cygnus Technologies is available to apply its proven technologies to develop such antibodies and assays on custom basis.

Principle of the Procedure

The *E. coli* assay is a two-site immunoenzymetric assay. Samples containing *E. coli* HCPs are reacted with a horseradish peroxidase (HRP) enzyme labeled anti-*E. coli* antibody simultaneously in microtiter strips coated with an affinity purified capture anti-*E. coli* antibody. The immunological reactions result in the formation of a sandwich complex of solid phase antibody-HCP-enzyme labeled antibody. The microtiter strips are washed to remove any unbound reactants. The substrate, tetramethylbenzidine (TMB) is then reacted. The amount of hydrolyzed substrate is read on a microtiter plate reader and is directly proportional to the concentration of *E. coli* HCPs present.

Reagents & Materials Provided

Component	Product #	
Anti- <i>E. coli</i> :HRP	F411	
Affinity purified antibody conjugated to HRP in a protein matrix with preservative. 1x12mL		
Anti-E. coli coated microtiter strips	F412*	
12x8 well strips in a bag with desiccant	1412	
E. coli HCP Standards	F413	
Solubilized E. coli HCPs in bovine serum		
albumin with preservative. Standards at 0, 1, 3,		
12, 40, and 100ng/mL. 1 mL/vial		
Stop Solution	F006	
0.5M sulfuric acid. 1x12mL		
TMB Substrate	F005	
3,3',5,5' Tetramethylbenzidine. 1x12mL		
Wash Concentrate (20X)	F004	
Tris buffered saline with preservative. 1x50mL		
*All components can be surphered concretely event # F112		

*All components can be purchased separately except # F412.

Materials & Equipment Required But Not Provided

- Microtiter plate reader spectrophotometer with dual wavelength capability at 450 & 650nm. (If your plate reader does not provide dual wavelength analysis you may read at just the 450nm wavelength.)
- Pipettors 25µL and 100µL
- Repeating or multichannel pipettor 100μL
- Microtiter plate rotator (400-600 rpm)

- Sample Diluent (recommended Cat # I028)
- Distilled water
- 1 liter wash bottle for diluted wash solution

Storage & Stability

- All reagents should be stored at 2°C to 8°C for stability until the expiration date printed on the kit.
- After prolonged storage, you may notice a salt precipitate and/or yellowing of the wash concentrate. These changes will not impact assay performance. To dissolve the precipitate, mix the wash concentrate thoroughly and dilute as directed in the 'Preparation of Reagents' section. Reconstituted wash solution is stable until the expiration date of the kit.

Precautions

- For Research or Manufacturing use only.
- Stop reagent is 0.5M H₂SO₄. Avoid contact with eyes, skin, and clothing.
- This kit should only be used by qualified technicians.

Preparation of Reagents

- Bring all reagents to room temperature.
- Dilute wash concentrate to 1 liter in distilled water, label with kit lot and expiration date, and store at 4°C.

Procedural Notes

1. Complete washing of the plates to remove excess unreacted reagents is essential to good assay reproducibility and sensitivity. We advise against the use of automated or other manually operated vacuum aspiration devices for washing plates as these may result in lower specific absorbances, higher non-specific absorbance, and more variable precision. The manual wash procedure described below generally provides lower backgrounds, higher specific absorbance, and better precision. If duplicate CVs are poor, or if the absorbance of the '0' standard is greater than 0.200, evaluate plate washing procedure for proper performance.

2. High Dose Hook Effect or poor dilutional linearity may be observed in samples with very high concentrations of HCP. High Dose Hook Effect is due to insufficient excess of antibody for very high concentrations of HCPs present in samples upstream in the purification process. Samples greater than 200µg/mL may give absorbances less than the 100ng/mL standard. It is also possible for samples to have certain HCPs in concentrations exceeding the amount of antibody for that particular HCP. In such cases the absorbance of the sample at all dilutions may be lower than the highest standard in the kit, however these samples will fail to show acceptable dilutional linearity/parallelism as evidenced by an apparent increase in dilution corrected HCP concentration with increasing dilution. High Dose Hook and poor dilutional linearity are most likely to be encountered from samples early in the purification process. If a hook effect is possible, samples should also be assayed diluted. If the HCP concentration of the undiluted sample is less than the diluted sample this may be indicative of the hook effect. Such samples should be diluted at least to the minimum required dilutions (MRDs) as established by your qualification studies using your actual final and in-process drug samples. The MRD is the first dilution at which all subsequent dilutions vield the same HCP value within the statistical limits of assav precision. The HCP value to be reported for such samples is the dilution corrected value at or greater than the established MRD. The diluent used should be compatible with accurate recovery. The preferred diluent is our Cat# I028 available in 100mL, 500mL, or 1 liter bottles. This is the same material used to prepare the kit standards. As the sample is diluted in 1028, its matrix begins to approach that of the standards, thus reducing any inaccuracies caused by dilutional artifacts. Other prospective diluents must be tested for non-specific binding and recovery by using them to dilute the 100ng/mL standard, as described in the "Limitations" section below

Limitations

- Before relying exclusively on this assay to detect host cell proteins, each laboratory should qualify that the kit antibodies and assay procedure yield acceptable specificity, accuracy, and precision. A suggested protocol for this qualification can be obtained from our Technical Services Department or our web site.
- The standards used in this assay are comprised of E. coli HCPs solubilized by mechanical disruption and detergent. 1D Western blot analysis of the antibodies used in this kit demonstrates that they recognize the majority of distinct PAGE separated bands seen using a sensitive protein staining method like silver stain or colloidal gold. Because the majority of HCPs will show sufficient antigenic conservation among all strains of E. coli this kit should be adequately reactive to HCPs from your strain. The antibodies used in this kit were generated against HCPs obtained predominately from DH5α and BL21 strains but they have been shown to react with the majority of HCP from all strains of E. coli tested. However, there can be no guarantee that this assay will detect all proteins or protein fragments from your process. If you desire a much more sensitive and specific method than western blot to detect the reactivity of the antibodies in this kit to your individual HCPs Cvanus is pleased to perform AAE as a service to

provide coverage information of the antibodies to the HCPs in your process samples.

- Certain sample matrices may interfere in this assay. The standards used in this kit attempt to simulate typical sample protein and matrices. However the potential exists that the product itself or other components in the sample matrix may result in either positive or negative interference in this assay. High or low pH, detergents, urea, high salt concentrations, and organic solvents are some of the known interference factors. It is advised to test all sample matrices for interference by diluting the 100ng/mL standard, 1 part to 4 parts of the matrix containing no or very low HCP impurities. This diluted standard when assaved as an unknown, should give an added HCP value in the range of 15 to 25 ng/mL. Consult Cygnus Technologies Technical Service Department for advice on how to quantitate the assav in problematic matrices.
- Avoid the assay of samples containing sodium azide (NaN₃) which will destroy the HRP activity of the conjugate and could result in the underestimation of HCP levels.

Quality Control

- Precision on duplicate samples should yield average % coefficients of variation of less than 10% for samples in the range of 3-100ng/mL. CVs for samples less than 3 ng/mL may be greater than 10%.
- It is recommended that each laboratory assay appropriate quality control samples in each run to ensure that all reagents and procedures are correct.

Calculation of Results

The standards may be used to construct a standard curve with values reported in ng/mL "total immunoreactive HCP equivalents". This data reduction may be performed through computer methods using curve-fitting routines such as point-to-point, cubic spline, or 4 parameter logistic fit. **Do not use linear regression analysis to interpolate values for samples as this may lead to significant inaccuracies!** Data may also be manually reduced by plotting the absorbance values of the standard on the y-axis versus concentration on the x-axis and drawing a smooth point-to-point line. Absorbances of samples are then interpolated from this standard curve.

Assay Protocol

 The assay is very robust such that assay variables like incubation times, sample size, and other sequential incubation schemes can be altered to manipulate assay performance for more sensitivity, increased upper analytical range, or reduced sample matrix interference. Before modifying the protocol from what is recommended, you are advised to contact Technical Services for input on the best way to achieve your desired goals.

- The protocol specifies use of an approved orbital microtiter plate shaker for the immunological steps. These can be purchased from most laboratory supply companies. If you do not have such a device, it is possible to incubate the plate without shaking however it will be necessary to extend the immunological incubation step in the plate by about one hour in order to achieve comparable results to the shaking protocol. Do not shake during the 30 minute substrate incubation step, as this may result in higher backgrounds and worse precision.
- Bring all reagents to room temperature. Set-up plate spectrophotometer to read dual wavelength at 450nm for the test wavelength and ~650nm for the reference.
- Thorough washing is essential to proper performance of this assay. Automated plate washing systems or other vacuum aspiration devices are not recommended. The manual method described in the assay protocol is preferred for best precision, sensitivity and accuracy. A more detailed discussion of this procedure can be obtained from our Technical Services Department or on our web site. In addition, a video demonstration of proper plate washing technique is available in the 'Technical Help' section of our web site.
- All standards, controls, and samples should be assayed at least in duplicate.
- Maintain a repetitive timing sequence from well to well for all assay steps to ensure that all incubation times are the same for each well.
- Make a work list for each assay to identify the location of each standard, control, and sample.
- It is recommended that your laboratory assay appropriate quality control samples in each run to ensure that all reagents and procedures are correct. You are strongly urged to make controls in your typical sample matrix using HCPs derived from your cell line. These controls can be aliquoted into single use vials and stored frozen for long-term stability.
- If the substrate has a distinct blue color prior to assay it may have been polluted. If the absorbance of 100μL of substrate plus 100μL of stop against a water blank is greater than 0.1 it may be necessary to obtain new substrate or the sensitivity of the assay may be compromised.
- Strips should be read within 30 minutes after adding stop solution since color will fade over time.

Assay Protocol

1. Pipette $25 \mu L$ of standards, controls and samples into wells indicated on work list.

2. Pipette 100µL of anti-*E. coli*:HRP (#F411) into each well.

3. Cover & incubate on orbital shaker at 400-600 rpm for 90 minutes at room temperature, $24^{\circ}C \pm 4^{\circ}C$.

4. Dump contents of wells into waste. Blot and gently but firmly tap over absorbent paper to remove most of the residual liquid. Overly aggressive banging of the plate or use of vacuum aspiration devices in an attempt to remove all residual liquid is not necessary and may cause variable dissociation of antibody bound material resulting in lower ODs and worse precision. Fill wells generously to overflowing with diluted wash solution using a squirt bottle or by pipetting in ~350µL. Dump and tap again. Repeat for a total of 4 washes. Wipe off any liquid from the bottom outside of the microtiter wells as any residue can interfere in the reading step. Do not allow wash solution to remain in wells for longer than a few seconds. Do not allow wells to dry before adding substrate.

5. Pipette 100µL of TMB substrate (#F005).

6. Incubate at room temperature for 30 minutes. DO NOT SHAKE.

7. Pipette 100µL of Stop Solution (#F006).

8. Read absorbance at 450/650nm.

Well #	Contents	Abs. at 450- 650nm	Mean Abs.
A1	Zero Std	0.066	0.067
B1	Zero Std	0.068	
C1	1ng/mL	0.099	0.099
D1	1ng/mL	0.098	
E1	3ng/mL	0.155	0.154
F1	3ng/mL	0.153	
G1	12ng/mL	0.396	0.205
H1	12ng/mL	0.393	0.395
A2	40ng/mL	1.190	1.185
B2	40ng/mL	1.179	
C2	100ng/mL	2.481	2.493
D2	100ng/mL	2.505	

Example Data

Performance Characteristics

Cygnus Technologies has qualified this assay by conventional criteria as indicated below. A copy of this qualification report can be obtained on our web site or by request. This gualification is generic in nature and is intended to supplement but not replace certain user and product specific qualification and qualification that should be performed by each laboratory. At a minimum each laboratory is urged to perform a spike and recovery study in their sample types. In addition, any of your samples types containing process derived HCPs within or above the analytical range of this assay should be evaluated for dilutional linearity to ensure that the assay is accurate and has sufficient antibody excess for your particular HCPs. Each laboratory and technician should also demonstrate competency in the assay by performing a precision study similar to that described below. A more detailed discussion of recommended user qualification protocols can be obtained by contacting our Technical Services Department or at our web site.

Sensitivity

The lower limit of detection (LOD) is defined as that concentration corresponding to a signal two standard deviations above the mean of the zero standard. LOD is ~0.2 ng/mL.

The lower limit of quantitation (LOQ) is defined as the lowest concentration, where concentration coefficients of variation (CVs) are less than 20%. The LOQ is ~1 ng/mL.

Specificity/Cross-Reactivity

1D Western blot and ELISA analysis against several strains of E. coli (DH5a, BL21, JM109, TOP10F, K12, & MC1061) indicate that most of the proteins are conserved among all strains. Thus, this assay should be useful for detecting HCPs from other E. coli cell lines. Each end user must qualify that this kit is adequately reactive and specific for their samples. Western blot, both 1 & 2 dimensional, is highly orthogonal to ELISA and to non-specific protein staining methods such as silver stain or colloidal gold. As such, the lack of identity between silver stain and western blot does not necessarily mean there is no antibody to that protein or that the ELISA will not detect that protein. If you desire a much more sensitive and specific method than Western blot to detect the reactivity of the antibodies in this kit to your individual HCPs Cygnus is pleased to perform AAE as a service to provide coverage information of the antibodies to the HCPs in your process samples. This method has been shown to be at least 100 fold more sensitive than Western blots in detecting antibody reactivity to individual HCPs. The same antibody as is used for both capture and HRP label can be purchased separately.

Cross reactivity to non-HCP components has not been extensively investigated with this kit. You should

evaluate components in your samples for positive interferences such as cross reactivity and non-specific binding. Negative interference studies are described below.

Precision

Both intra (n=20 replicates) and inter-assay (n=5 assays) precision were determined on 3 pools with low (~3ng/mL), medium (~12ng/mL), and high concentrations (~40ng/mL). The % CV is the standard deviation divided by the mean and multiplied by 100.

Pool	Intra assay CV	Inter assay CV
Low	3.3%	7.7%
Medium	1.9%	5.5%
High	2.7%	6.5%

Recovery/Interference Studies

Various buffer matrices were evaluated by adding known amounts of E. coli HCP preparation used to make the standards in this kit. Because this assay is designed to minimize matrix interference most of these buffers vielded acceptable recovery defined as between 80-120%. The standards used in this kit contain 8mg/mL of bovine serum albumin intended to simulate non-specific protein effects of most sample proteins or pDNA products. However, very high concentrations of some products (often in the 2-5 mg/mL range) may interfere in the accurate measurement of HCPs. In general, extremes in pH (less than 5.0 and greater than 8.5), high salt concentration, high polysaccharide concentrations, urea, organic solvents, and most detergents can cause under-recovery. Each user should qualify that their sample matrices yield accurate recovery. Such an experiment can be performed, by diluting the 100ng/mL standard provided with this kit, into the sample matrix in question as described in the "Limitations" section. Cvanus offers a more concentrated form of the HCP used to prepare the kits standards for your spike recovery and preparation of analyte controls.

Hook Capacity

Increasing concentrations of HCPs greater than 100 ng/mL were assayed as unknowns. The hook capacity, defined as that concentration yielding an absorbance reading less than the 100 ng/mL standard was greater than 200 μ g/mL.

Ordering Information/ Customer Service

Cygnus Technologies also offers kits for the extraction and detection of *E. coli* Host Cell DNA. The following kits are available:

Residual Host Cell DNA extraction:

Cat # D100W, DNA Extraction Kit in 96 deep well plate Cat # D100T, DNA Extraction Kit in microfuge tubes

 Extraction and PCR amplification of *E. coli* Host Cell DNA for use with user supplied master mix:

Cat # D415W, DNA Extraction Kit in 96 deep well plate Cat # D415T, DNA Extraction Kit in microfuge tubes

• Residual *E. coli* Host Cell DNA extraction and detection using PicoGreen® dye:

Cat # D410W, DNA Extraction Kit in 96 deep well plate Cat # D410T, DNA Extraction Kit in microfuge tubes

To place an order or to obtain additional product information contact *Cygnus Technologies*:

www.cygnustechnologies.com Cygnus Technologies, LLC 4332 Southport Supply Rd. SE Southport, NC 28461 USA Tel: 910-454-9442 Fax: 910-454-9443 Email: techsupport@cygnustechnologies.com